



## **MLGA (MGA) in Feed**

**January 2010 R-Biopharm Inc**

**R-Biopharm AG**  
**Landwehrstraße 54**  
**64293 Darmstadt**

**Telefon: 0 61 51 / 81 02-0**  
**Telefax: 0 61 51 / 81 02-40**  
**info@r-biopharm.de**  
**www.r-biopharm.com**

## MLGA

### RIDASCREEN MELENGESTROL ACETATE (MLGA)

- ◆ Competitive Enzyme Immunoassay
- ◆ Sample Extraction: Homogenize, Dilute, Centrifuge, Solid-phase Cleanup- (Approx 2.5 hours per sample.)
- ◆ Incubation time 2.5 hours
- ◆ Cross-Reactivity: Melengestrol Acetate 100%, Megestrol Acetate 10%, Medroxyprogesterone Acetate 6.6%,

## MLGA

### Overview

- **Synthetic gestagens:** Used for estrus inhibition or synchronization.
- **Melengestrol acetate** belongs to the most active synthetic gestagens.
- **Oral bioactivity:** 10 to 100 times higher than gestagens chlormadinone acetate (CMA) or medroxyprogesterone acetate.
- **MLGA** is a licensed growth promoting feed additive for heifers in the USA and Canada.
- **Admitted dose** is 0.5 mg per day and head, given as a feed premix.
- **Mode of anabolic action** is unclear, but it stimulates the ovarian synthesis of the endogenous anabolic steroid estradiol which may have androgenic side effects.
- **Parenterally**, hormonal activity of MLGA is still 125 times higher than that of progesterone.
- **Strongly lipophilic**, MLGA is accumulated in fat 200x higher than in blood plasma.
- **European Union:** Use of sexual hormones for growth promoting purposes is generally forbidden since 1988, as well as the import of meat from hormone treated cattle.

## MLGA

### Extraction for Feed and Premixes

- Homogenize approximately 25 g of feed sample.
- Transfer 1.0 g of the homogenized sample into a glass tube.
- Add 7.0 mL of 80 % methanol in distilled water.
- Mix for 30 sec using a vortex.
- Mix for another 10 min using a shaker.
- Add 4.0 mL of distilled water and mix again for 30 sec using a vortex.
- Centrifuge sample for 6 minutes at 4000 rpm.
- Transfer 5.5 mL of the clean extract onto an activated C18 cartridge.
- Conduct the solid phase clean-up procedure

**MLGA**

## **Solid Phase Clean-Up**

- Activate solid phase cartridges (C18, R-Biopharm Cat. No. R2002)
  - wash the cartridges subsequently with 5 mL MeOH 2x and with 5 ml 50 % MeOH in distilled water two time.
- Transfer 5.5 mL of feed sample extract onto activated solid phase cartridge.
- Wash the cartridges using 7.0 mL 50 % methanol in distilled water.
- Dry the cartridges under vacuum.
- Elute using 1 ml 80 % methanol in distilled water with a speed of 1 mL per min.
- Dilute extract 1+1 with distilled water.
- Use 20 µL per well in the assay.

Depending on the MGA concentration in the sample, further dilutions may be necessary. In this case use methanol/distilled water (40/60 v/v) for all further dilutions (i.e. always maintain 40% methanol)

## MLGA

### Test Procedure

- Insert a sufficient number of wells into the microwell holder for all standards and samples to be run in duplicate.
  - Add 20  $\mu$ L of each standard solution or prepared sample to separate duplicate wells.
  - Add 50  $\mu$ L of diluted enzyme conjugate to each well.
  - Add 50  $\mu$ L of the diluted anti-MLGA antibody solution to each well.
- Mix and incubate 2 h at room temperature.
- Pour the liquid out of the wells and rinse 3x with wash solution.
  - Add 50  $\mu$ L of substrate and 50  $\mu$ L of chromogen to each well. Mix and incubate for 30 min at room temperature in the dark.
  - Add 100  $\mu$ L of the stop solution to each well.
  - Mix and measure the absorbance at 450 nm.

**MLGA**

**Results:**

Initial studies were performed using feed and premix samples provided by SDSU.

Subsequent data was created with a commercially available feed, with and without addition of a premix. LOD was calculated using this feed only.

## R-Biopharm Group

Argentina • Austria • Brazil • China  
France • Germany • Italy • Canada  
Netherlands • Switzerland • UK • USA



MLGA

MGA	Claim (ppb)	Spike added (ppb)	Result	Dilutions	RPD & spike recovery
09F-04167A	2200		3743	(1:1250)	4.1
09F-04167B	2200		3594	(1:1250)	
09F-04167C (spike)	2200	2000	5261	(1:1250)	94%
09F-04167D (spike)	2200	2000	5358	(1:1250)	
09F-03896A	882				
09F-03896B (spike)	882	1000			
08S-20786A	881		824	(1:625)	14.6
08S-20786B	881		712	(1:625)	
08S-20786C (spike)	881	1000	1540	(1:625)	91%
08S-20786D (spike)	881	1000	1661	(1:625)	
Steak-Maker (40-28) A	880		364	(1:625)	8.3
Steak-Maker (40-28) B	880		335	(1:625)	
Steak-Maker (40-28) C (spike)	880	1000	965	(1:625)	73%
Steak-Maker (40-28) D (spike)	880	1000	1010	(1:625)	
11 ppm Spiking Solution	11000		11190	(1:6250)	10.7
12 ppm Spiking Solution	11001		10050	(1:25000)	
Premix A	220000		296800	(1:125000)	2.5
Premix B	222000		289600	(1:125000)	
Premix C Spike	220000	110000	376000	(1:178600)	97%
Premix D Spike	222000	110000	403857	(1:178600)	



**MLGA**

## **SDSU Summary:**

- **RPD (Range):  
2.5-14.6**
- **RPD (Mean):  
8.0**
- **Percent Recovery (Range):  
73%-97%**
- **Percent Recovery (Mean):  
89%**

## R-Biopharm Group

Argentina • Austria • Brazil • China  
France • Germany • Italy • Canada  
Netherlands • Switzerland • UK • USA



### MLGA

Sample	Claim	Dilution	Replicates	Results	
With MLGA	880 ppb	1:625	20	Mean	845.35
				SD	47.27
				CV	5.59
No MLGA	0 ppb	1:4	20	Mean	0.62
				SD	0.03
				CV	5.52
				LOD	0.72
Premix	220000 ppb	1:125000	10	Mean	228730.00
				SD	3369.82
				CV	1.47

**MLGA**

**References:**

1. Daxenberger, A. et al.; Vet. Quart. 21, 154-158 (1999):

Detection of Melengestrol acetate residues in plasma and edible tissues of Heifers

2. SCREENING OF MELENGESTROL ACETATE (MGA) IN FAT AND MUSCLE BY ENZYME IMMUNOASSAY (EIA)

Melanie Hageleit, Andreas Daxenberger and Heinrich H.D. Meyer

Institut für Physiologie, D - 85350 Freising-Weihenstephan, Germany

## R-Biopharm Group

Argentina • Austria • Brazil • China  
France • Germany • Italy • Canada  
Netherlands • Switzerland • UK • USA



MLGA

# Thank you for your attention!



[www.r-biopharm.com](http://www.r-biopharm.com)